2017-12-14 / HS

Produktmanual AC 213500

English
manual
overleaf

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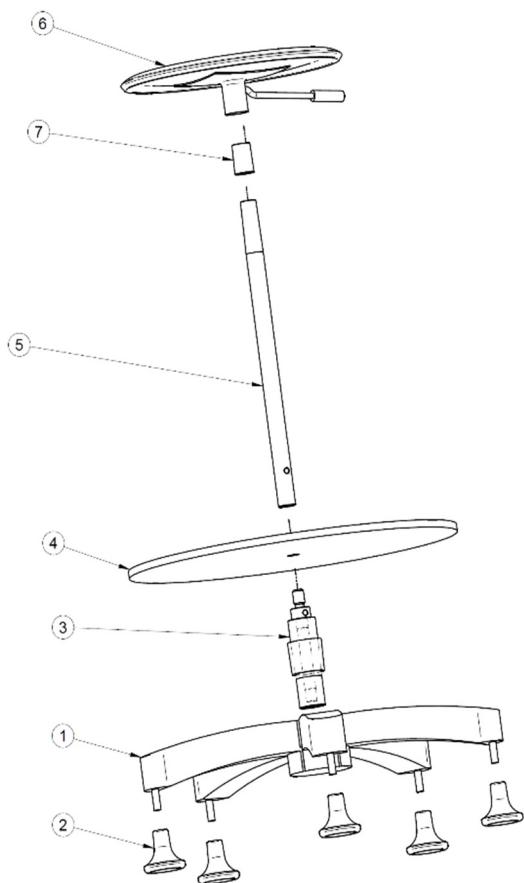
The rotating platform is used, e.g. with bicycle wheel gyro 213600, to demonstrate conservation of angular momentum.

Assembly

- I. Mount the feet (2) under the base (1)
- II. Insert the cylinder (3) into the base with the threading facing upwards
- III. Place the foot rest (4) with the threading of the cylinder (3) through the centre hole.
- IV. Screw the rod (5) onto the cylinder (3)
- V. Insert the short grey plastic tube (7) below the seat
- VI. Place the seat (with plastic tube) on top of the rod

Applications – I

Select preferably a small and light person. (However, the person must have arms that are long and strong enough to handle the gyro as described below.)

The test person sits down on the stool on the rotating platform and holds the bicycle wheels in stretched arms. Start for example with the shaft vertically with the right hand at the top.

The test person grips the handles firmly while a helper puts the gyro in fast motion.

Still with stretched arms, the test person rotates the gyro's pivot axis until the left hand comes to the top. The platform and the test person will now rotate slowly in the original direction of rotation of the gyro.

Applications – II

The test person sits down on the stool on the rotating platform and holds two dumbbells in stretched arms. A helper makes the test person rotate **slowly**.

The test person moves the dumbbells together, close to the chest. The rotational speed increases significantly.

Caution!

In case of careless use,
the test person may fall off.

In the case of careless use of the bicycle wheel gyro, it can scrape against the forearms.

Frederiksen Scientific cannot be held liable for damage caused by careless use of the equipment.

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© Frederiksen Scientific A/S