280050_AC

21.11.16

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21.11.16

Description

This experiment lamp is a multi-purpose source of white light based on a halogen bulb. The lamp is provided with a cooling fan to keep the temperature of the housing below 50 °C. The filament position can be adjusted both longitudinally and laterally. The entire lamp can be angled vertically. An optional holder for apertures can be mounted in front of the lens. The lamp housing is fitted with a Ø10mm support rod and 2 safety sockets for connection to the power supply.

Required accessories

- 0006.00 Stand base.
- 3610.55/3618.60/3630.00 Power supply.
- 2 test leads

Additional accessories

- 2800.55 Aperture holder for experiment lamp.

Operating instructions

Connect the lamp to a power supply with two test leads. The voltage must not exceed 12 V AC/DC!

Adjusting the lamp: Loosen the screw below the adjustment handle. The socket for the bulb can now be adjusted in the longitudinal direction by pulling / pushing the handle. Likewise, the socket is adjusted laterally inside the house by rotating the handle. When the filament is in the desired position it can be fixed by tightening the thumbscrew. The lamp housing can be tilted vertically by means of the fixing screw below the lamp housing. Replacing the halogen bulb: Remove the connection to the power supply. Remove the 4 thumbscrews at the top of the front and back plates. Lift off the lid. You now have access to change the lamp.

IMPORTANT:

When inserting the halogen lamp, the socket should be pushed as far forward toward the lens as possible. This avoids the risk of bending the metal plate on which the socket is mounted which would hinder the free movement of the socket. Install the cover again before use. Note that the vents should be at the front of the lid.

Spare parts

Halogen lamp 12 V / 50 W (4270.10)

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© A/S Søren Frederiksen, Ølgod