## Geiger-Müllererrør på stang med BNC-stik nr. 5125.15

06.07.12

AC 5125.15

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06.07.12

AC 5125.15

The GM-tube no. 5125.15 is well suited for measuring alpha, beta and gamma radiation.

For mounting the GM tube is equipped with a steel rod 10 mm in diameter.

The GM-tube no. 5125.15 comes complete with cable mounted with a BNC plug for connection to GM counters with BNC socket.

The tube's window is made of a very thin and delicate material which is easily destroyed. For this reason the tube is delivered with a protective cap. It is recommended always to use this cap for protection when the tube is not in use. The cap is provided with a venting hole to avoid vaccum when removing it or the opposit when mounting it again. I.e. it is of the utmost importance that this venting hole is uncovered when mounting or removing the cap. For certain applications removing the cap is quite unnecessary e.g. gamma rays will penetrate the cap without any problems.

**Note:** Although the GM tube itself is reasonably well protected in the plastic housing, it is still a fragile component. Do not drop, hit or squeeze the house. Use the included steel rod for mounting.t

Please notice that the GM-tube has no built-in anode resistor. Our GM counters are provided with a built-in anode resistor.

### Technical specifications:

The tube is sensitive to alpha, beta and gamma radiation.

The length of the tube is 32 mm.

The window is retracted 4,5 mm from the front of the housing.

Plateau threshold voltage ( $V_{b1}$ ): 450 V

Plateau length ( $V_{b2} - V_{b1}$ ): 150 V

Recommended supply voltage: 500 V

### Window specifications:

Thickness: 1.5 til 2.0 mg/cm<sup>2</sup>

Usefull diameter: 9.1 mm

Material: Mica

Gas filling: Neon and Halogen

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© A/S Søren Frederiksen, Ølgod